


HPLC Method Limited Verification for TGF-001 (non-GMP)


Project	Pyronaridine_INV-054926
Compound	TGF-001
Purpose	Method Verification
Category	Methods
Substance Type	Drug Substance
Report ID	INV_054926_HPLC_V4 Version 1.0

Authors

Cai Alain (Alain)

Approvers

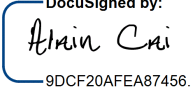By:  9DCF20AFEA87456... Title: AD Manager Date: 1/21/2024

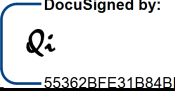By:  55362BFE31B84BD... Title: Head of Technical Development Date: 1/21/2024

Distribution

Zhou Irvin, Pyronaridine (INV-054926) project team, Approvers

Table of contents


1.	Objective	3
2.	Summary and conclusion	3
3.	Results	6
3.1.	Autosampler stability	6
3.2.	Specificity/ selectivity	7
3.3.	LOD/ LOQ	9
3.4.	Linearity of TGF-001	13
3.5.	Linearity of DIA and DIN (RRF determination)	15
3.6.	Accuracy of TGF-001	17
3.7.	Repeatability	18
4.	Experimental	19
4.1.	Equipment	19
4.2.	Equipment and reagents	19
4.3.	Solutions	19
4.3.1.	Diluent	19
4.3.2.	Mobile phase	20
4.3.3.	Solutions	20

1. Objective

TGF-001 is analyzed by the newly developed HPLC method 'INV_054926_HPLC_M4' using a reversed phase C18 column, e.g.: Waters X-Bridge column. In this document, the limited verification of the new HPLC method shall be performed at non-GMP level.

Related reports:

INV_054926_HPLC_M4: HPLC Method Description for In-process control, Identity, Assay and Related Substances of TGF-001

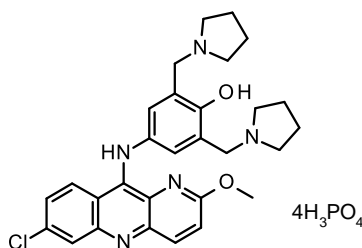

TGF-001

2. Summary and conclusion

The method verification has been performed at non-GMP level. The following analytical parameters have been assessed:

- Autosampler stability
- Specificity/ selectivity
- LOD
- LOQ
- Linearity of TGF-001
- Linearity of DIA and DIN (RRF determination)
- Accuracy of TGF-001
- Repeatability

The peak purity showed that unknown peaks at RRT 0.85/ 0.87 should be coeluted peaks. The obtained results are summarized in Table 1.

Table 1 non-GMP verification of the new developed HPLC method for TGF-001

Parameter	Acceptance criteria	Results	Conformity	Remark
Autosampler stability	For TGF-001: % Δ \leq 2% For impurities (report all \geq LOQ): between LOQ and \leq 0.1%: % Δ \leq 40%, > 0.1% and \leq 0.5%: % Δ \leq 20%, > 0.5%: % Δ \leq 15% Report limit time for autosampler stability	Stable at least 24 hours for assay Sample needs freshly prepared for impurity quantification	Pass	See section 3.1
Specificity/ selectivity	No interference between the blank peaks and the components of interest in blanks. All impurity peaks should be separated with TGF-001 peak	No interference peaks in blanks. All impurity peaks are separated with TGF-001 peak.	Pass	Unknown peaks at RRT 0.85/0.87 should be coeluted peaks. See section 3.2
LOD	S/N (0.03% TGF-001) \geq 3 S/N (0.03% DIA) \geq 3 S/N (0.03% DIN) \geq 3	S/N (0.03% TGF-001) = 38.2 S/N (0.03% DIA) = 37.6 S/N (0.03% DIN) = 60.4	Pass	See section 3.3
LOQ	S/N (0.05% TGF-001) \geq 10 S/N (0.05% DIA) \geq 10 S/N (0.05% DIN) \geq 10 %RSD (0.05% TGF-001 peak area) \leq 30 % %RSD (0.05% DIA peak area) \leq 30 % %RSD (0.05% DIN peak area) \leq 30 %	S/N (0.05% TGF-001) = 60.5 S/N (0.05% DIA) = 73.0 S/N (0.05% DIN) = 83.2 %RSD (0.05% TGF-001) = 1.3% %RSD (0.05% DIA) = 0.9% %RSD (0.05% DIN) = 0.2%	Pass	See section 3.3
Linearity of TGF-001	a) LOQ to 140%: $R^2 \geq 0.995$, y intercept \leq 5% of the nominal concentration, Report slope, show figure. Report R^2 , y intercept, slope, figure for: b) LOQ to 5% c) 50% to 140%	LOQ to 140%: $R^2 = 0.9999$ y intercept = 0.37% Slope = 52689 LOQ to 5%: $R^2 = 0.9998$ y intercept = 0.05% Slope = 51576 50% to 140%:	Pass	See section 3.4

		$R^2 = 0.9998$ y intercept = 2.19% Slope = 53579		
Linearity of DIA and DIN (RRF determination)	LOQ to 0.5%: $R^2 \geq 0.995$, Report slope, show figure. Calculate RRF value.	DIA: $R^2 = 0.9995$ Slope = 45379 RRF = 1.16 DIN: $R^2 = 0.9993$ Slope = 58071 RRF = 0.91	Pass	See section 3.5
Accuracy of TGF-001	%RSD $\leq 2\%$ for each concentration level Single recovery: 95% - 105% Mean recovery for each concentration level: 98% - 102% (Recovery is calculated as determined concentration vs. nominal concentration)	%RSD (80%) = 0.6% %RSD (100%) = 1.7% %RSD (120%) = 2.0% Recovery (80% prep. 1) = 99% Recovery (80% prep. 2) = 99% Recovery (80% prep. 3) = 98% Recovery (100% prep. 1) = 100% Recovery (100% prep. 2) = 100% Recovery (100% prep. 3) = 97% Recovery (120% prep. 1) = 101% Recovery (120% prep. 2) = 98% Recovery (120% prep. 3) = 98% Mean recovery (80%) = 98% Mean recovery (100%) = 99% Mean recovery (120%) = 99%	Pass	See section 3.6
Repeatability	For TGF-001: %RSD (100% TGF-001, n=6) $\leq 2\%$ For DIA: %RSD (0.14% DIA, n=6) $\leq 15\%$ For DIN: %RSD (0.14% DIN, n=6) $\leq 15\%$	%RSD (100% TGF-001) = 0.1% %RSD (0.14% DIA in 100% TGF-001) = 0.6% %RSD (0.14% DIN in 100% TGF-001) = 0.9%	Pass	See section 3.7

3. Results

3.1. Autosampler stability

Autosampler stability was performed using TGF-001 batch 100144-200502-REF.

The sample solution was found to be stable at least 24 hours for assay purpose and instable for unknown impurity at RRT 0.82/ 0.85 which means fresh sample preparation is needed for impurity quantification.

Table 2 Results for autosampler stability of TGF-001

Time		0h			1h		2h		3h		4h		6h		9h		12h		18h		24h		Limit
RT, min	RRT	area%	area%	%Δ	area%	%Δ	area%	%Δ	area%	%Δ	area%	%Δ	area%	%Δ	area%	%Δ	area%	%Δ	area%	%Δ	area%	%Δ	%Δ
9.81	0.82	0.19%	0.24%	26%	0.26%	37%	0.28%	47%	0.29%	53%	0.31%	63%	0.32%	68%	0.33%	74%	0.34%	79%	0.35%	84%			≤ 20%
10.17	0.85	0.17%	0.15%	12%	0.14%	18%	0.13%	24%	0.13%	24%	0.12%	29%	0.12%	29%	0.11%	35%	0.12%	29%	0.11%	35%			≤ 20%
10.40	0.87	0.05%	0.05%	0%	0.04%	20%	0.04%	20%	0.04%	20%	0.04%	20%	0.04%	20%	0.03%	40%	0.03%	40%	0.03%	40%			≤ 40%
10.96	0.91	0.04%	0.04%	0%	0.04%	0%	0.04%	0%	0.04%	0%	0.04%	0%	0.05%	25%	0.05%	25%	0.05%	25%	0.05%	25%			-
11.24	0.94	0.05%	0.05%	0%	0.05%	0%	0.05%	0%	0.05%	0%	0.05%	0%	0.05%	0%	0.05%	0%	0.05%	0%	0.05%	0%			≤ 40%
11.33	0.95	0.05%	0.06%	20%	0.05%	0%	0.05%	0%	0.05%	0%	0.05%	0%	0.05%	0%	0.05%	0%	0.05%	0%	0.05%	0%			≤ 40%
11.99	1.00	98.25%	98.19%	0%	98.20%	0%	98.19%	0%	98.17%	0%	98.18%	0%	98.16%	0%	98.16%	0%	98.14%	0%	98.15%	0%			≤ 2%
12.80	1.07	0.34%	0.35%	3%	0.33%	3%	0.33%	3%	0.33%	3%	0.33%	3%	0.33%	3%	0.33%	3%	0.33%	3%	0.32%	6%			≤ 20%
16.77	1.40	0.08%	0.07%	13%	0.08%	0%	0.08%	0%	0.08%	0%	0.08%	0%	0.08%	0%	0.08%	0%	0.08%	0%	0.09%	13%			≤ 40%
23.46	1.96	0.09%	0.09%	0%	0.09%	0%	0.09%	0%	0.09%	0%	0.09%	0%	0.09%	0%	0.09%	0%	0.10%	11%	0.10%	11%			≤ 40%
24.83	2.07	0.52%	0.52%	0%	0.52%	0%	0.51%	2%	0.52%	0%	0.52%	0%	0.51%	2%	0.52%	0%	0.52%	0%	0.52%	0%			≤ 15%
29.92	2.50	0.20%	0.20%	0%	0.20%	0%	0.20%	0%	0.20%	0%	0.20%	0%	0.20%	0%	0.20%	0%	0.20%	0%	0.20%	0%			≤ 20%

$$\% \Delta = (\% \text{area}_{t=n} - \% \text{area}_{t=0}) / \% \text{area}_{t=0} * 100\%$$

3.2. Specificity/ selectivity

Specificity/ selectivity was performed using TGF-001 batch 100144-200502-REF and TGF-001 spiked with impurity DIA or DIN.

There is no interference between the blank peaks and the components of interest. All impurity peaks are separated with TGF-001 peak. The peak purity showed that unknown peaks at RRT 0.85/ 0.87 should be coeluted peaks.

Table 3 Results for specificity/ selectivity of TGF-001

TGF-001, batch		100144-200502-REF			
Name	RT, min	RRT	Area	% Area	peak purity
	9.806	0.82	16.1742	0.19%	995.4
	10.166	0.85	14.0179	0.17%	844.8
	10.404	0.87	4.5063	0.05%	618.1
	10.962	0.91	2.9933	0.04%	922.1
	11.240	0.94	3.9864	0.05%	988.4
	11.333	0.95	4.3823	0.05%	973.4
TGF-001	11.988	1.00	8336.9955	98.25%	1000.0
	12.803	1.07	28.4520	0.34%	997.0
	16.768	1.40	6.5751	0.08%	961.9
	23.464	1.96	7.4326	0.09%	989.1
	24.830	2.07	43.7429	0.52%	999.5
	29.915	2.50	16.5870	0.20%	999.3

Figure 1 Example HPLC chromatogram of blank, 278 nm

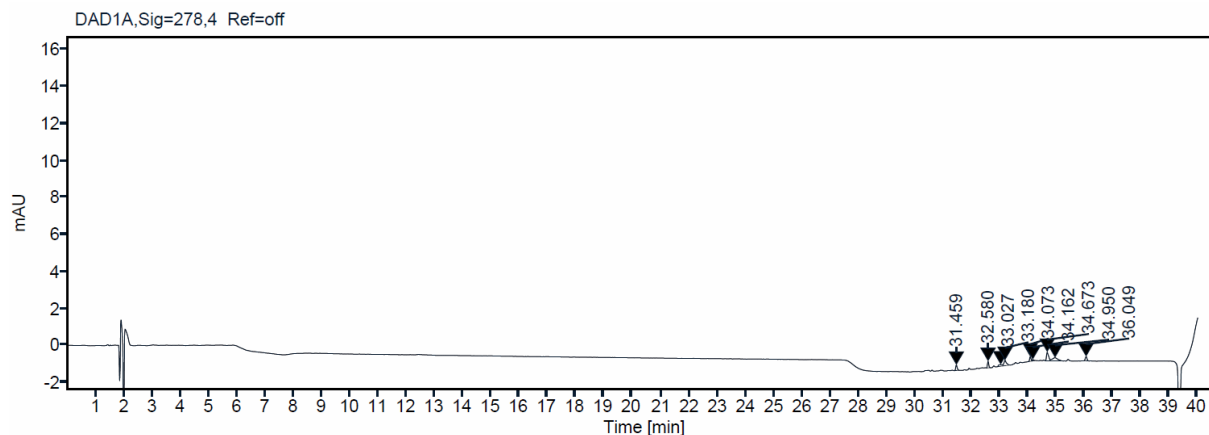
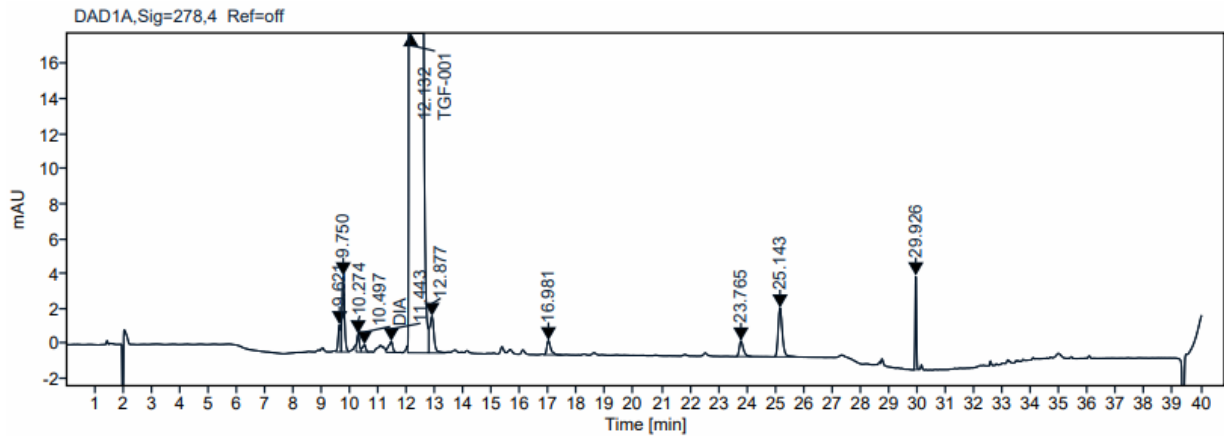
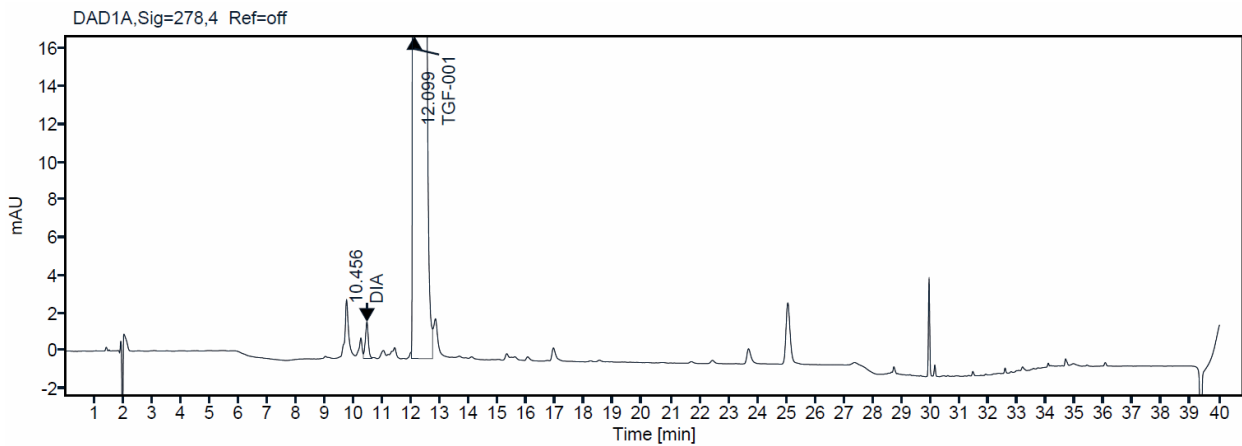
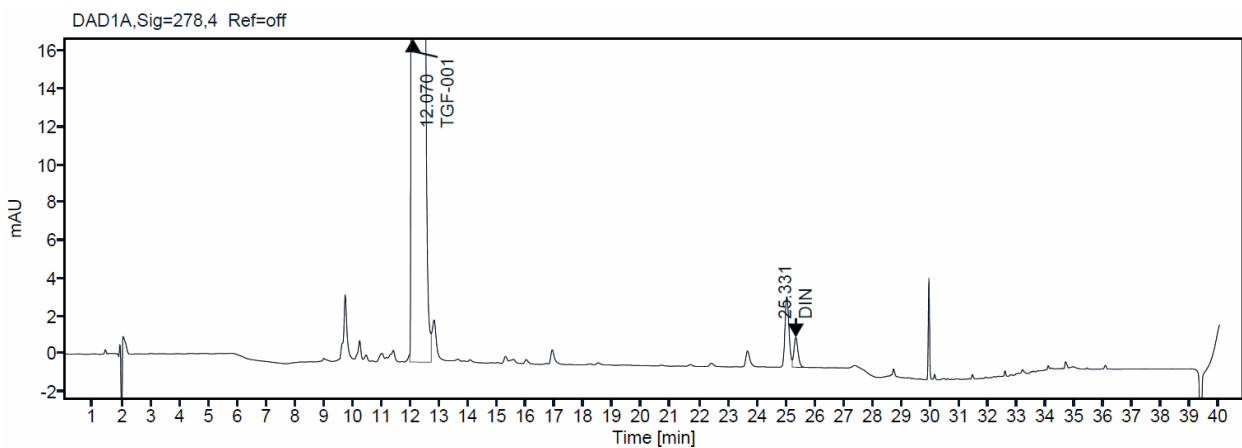


Figure 2 Example HPLC chromatogram of TGF-001 batch 100144-200502-REF, 278 nm**Figure 3** Example HPLC chromatogram of TGF-001 batch 100144-200502-REF spiked with impurity DIA batch PHTRACKD-394-REF, 278 nm**Figure 4** Example HPLC chromatogram of TGF-001 batch 100144-200502-REF spiked with impurity DIN batch PHTRACKD-389-REF, 278 nm

3.3. LOD/ LOQ

LOD and LOQ were performed using TGF-001 batch 100144-200502-REF, DIA batch PHTRACKD-394-REF and DIN batch PHTRACKD-389-REF.


The signal to noise of TGF-001 peak is calculated to be 38.2 in LOD solution (0.03%) and 60.5 in LOQ solution (0.05%) respectively. The %RSD of TGF-001 peak area in LOQ solutions (n=3) is found to be 1.3%.

The signal to noise of DIA peak is calculated to be 37.6 in LOD solution (0.03%) and 73.0 in LOQ solution (0.05%) respectively. The %RSD of DIA peak area in LOQ solutions (n=3) is found to be 0.9%. The signal to noise of DIN peak is calculated to be 60.4 in LOD solution (0.03%) and 83.2 in LOQ solution (0.05%) respectively. The %RSD of DIN peak area in LOQ solutions (n=3) is found to be 0.2%. All results meet the acceptance criteria.

Table 4 %RSD results for 3 injections of 0.05% solution

Reference solution 2	Area counts [pA]		
	0.05% TGF-001	0.05% DIA	0.05% DIN
Inj. 1	2.8677	2.3974	4.6605
Inj. 2	2.8049	2.3868	4.6678
Inj. 3	2.8068	2.4287	4.6801
Average	2.8265	2.4043	4.6695
STDEV	0.0357	0.0218	0.0099
%RSD	1.3%	0.9%	0.2%
Acceptance criteria	≤ 30%		


Figure 5 Example Chromatogram of signal to noise calculation for 0.03% TGF-001 solution (LOD).

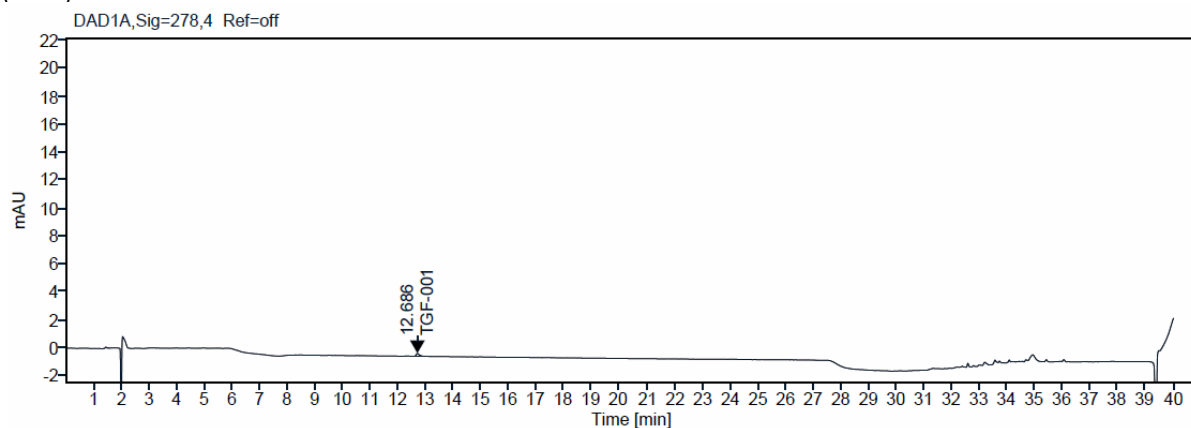

Signal: DAD1A, Sig=278,4 Ref=off

RT [min]	Type	Width [min]	Area	Height	Area%	Name	Peak Signal To Noise
12.686	MM m	0.3773	1.5783	0.1995	100.00	TGF-001	38.15009


Figure 6 Example chromatogram of signal to noise calculation for 0.05% TGF-001 solution (LOQ).

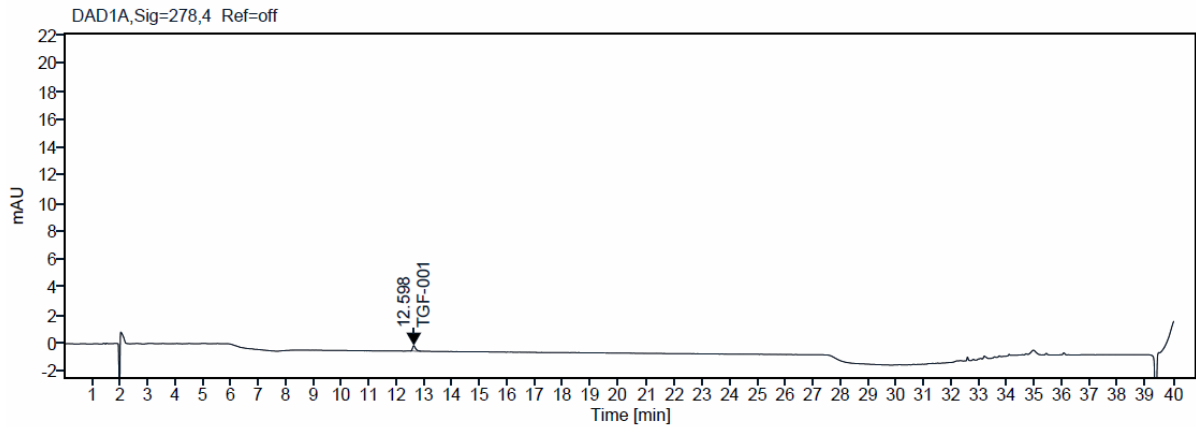

Signal: DAD1A,Sig=278,4 Ref=off


RT [min]	Type	Width [min]	Area	Height	Area% Name	Peak Signal To Noise
12.598	MM m	0.3746	2.8677	0.3789	100.00 TGF-001	60.546 25

Figure 7 Example Chromatogram of signal to noise calculation for 0.03% DIA solution (LOD).

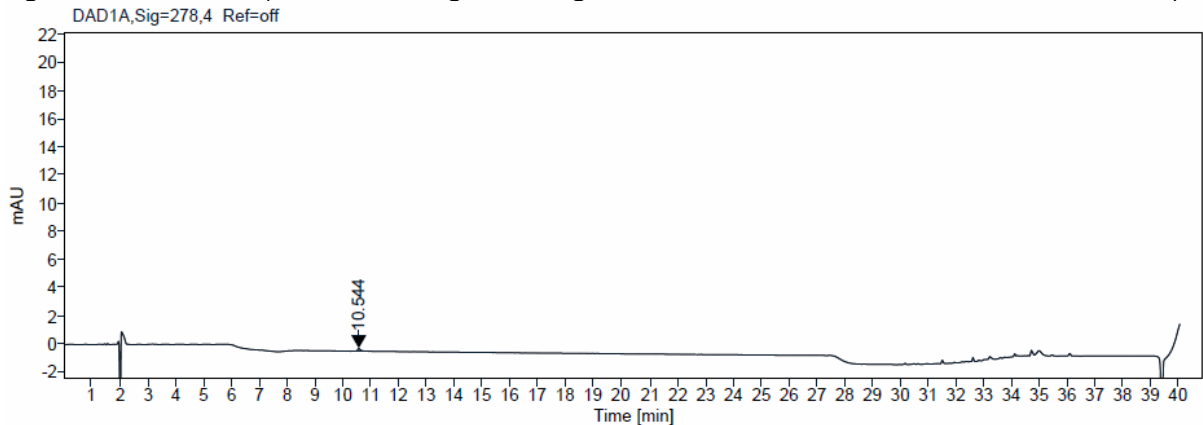


Signal: DAD1A,Sig=278,4 Ref=off

RT [min]	Type	Width [min]	Area	Height	Area% Name	Peak Signal To Noise
10.544	MM m	0.4546	1.4910	0.1968	100.00	37.566 24

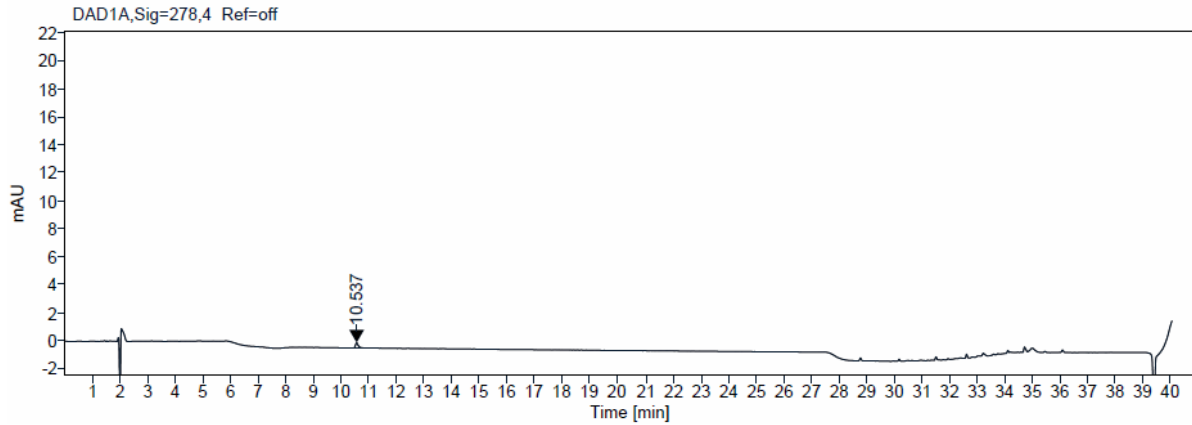Figure 8 Example Chromatogram of signal to noise calculation for 0.05% DIA solution (LOQ).


Signal: DAD1A,Sig=278,4 Ref=off

RT [min]	Type	Width [min]	Area	Height	Area%	Name	Peak Signal To Noise
10.537	MM m	0.3644	2.3974	0.3585	100.00		72.965 22

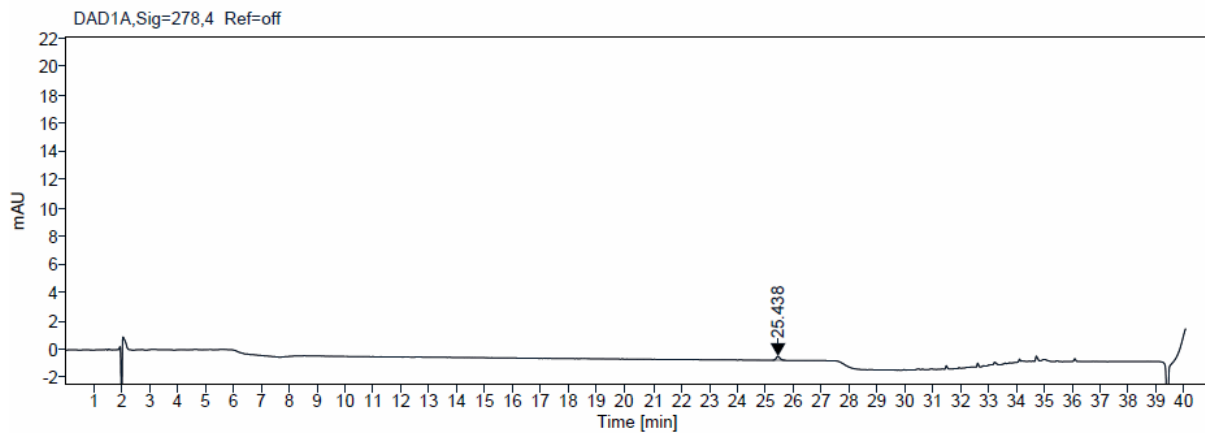Figure 9 Example Chromatogram of signal to noise calculation for 0.03% DIN solution (LOD).

Signal: DAD1A,Sig=278,4 Ref=off

RT [min]	Type	Width [min]	Area	Height	Area%	Name	Peak Signal To Noise
25.438	MM m	0.4408	2.5680	0.2710	100.00		60.431 99

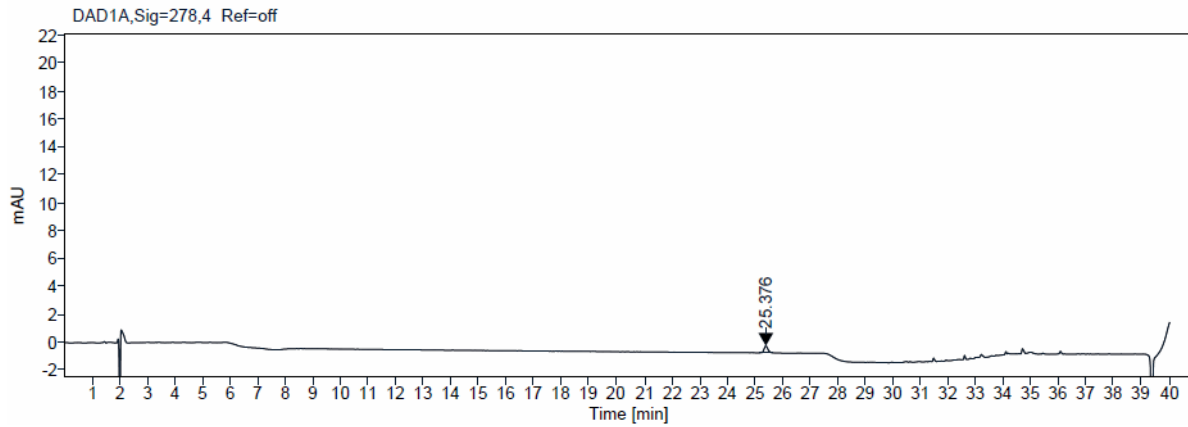Figure 10 Example Chromatogram of signal to noise calculation for 0.05% DIN solution (LOQ).

Signal: DAD1A,Sig=278,4 Ref=off

RT [min]	Type	Width [min]	Area	Height	Area%	Name	Peak Signal To Noise
25.376	MM m	0.4515	4.6605	0.4860	100.00		83.166 14

3.4. Linearity of TGF-001

Linearity was performed using TGF-001 batch 100144-200502-REF.

The correlation coefficient (R^2) was found to be 0.9999 for the range from LOQ to 140%, 0.9998 for the range from LOQ to 5% and 0.9998 for the range from 50% to 140%. The y intercept and slope are reported in Table 5. All results meet the acceptance criteria.

Table 5 Results for linearity of TGF-001

Items	Result	Acceptance criteria
LOQ to 140%		
R^2	0.9999	≥ 0.995
y intercept	0.37%	$\leq 5\%$ of the nominal concentration
Slope	52689	report
LOQ to 5%		
R^2	0.9998	report
y intercept	0.05%	report
Slope	51576	report
50% to 140%		
R^2	0.9998	report
y intercept	2.19%	report
Slope	53579	report

Table 6 Concentration and area counts of linearity for TGF-001

% of the conc.	Conc. (mg/mL)	Area counts [pA]
0.05%	0.00008	2.879
0.1%	0.00017	5.636
0.5%	0.00083	34.856
5%	0.00829	423.865
20%	0.03315	1682.543
50%	0.08288	4294.905
80%	0.13261	6852.764
100%	0.16576	8660.748
120%	0.19891	10518.338
140%	0.23207	12242.546

Figure 11 Linearity of TGF-001 from LOQ to 140%

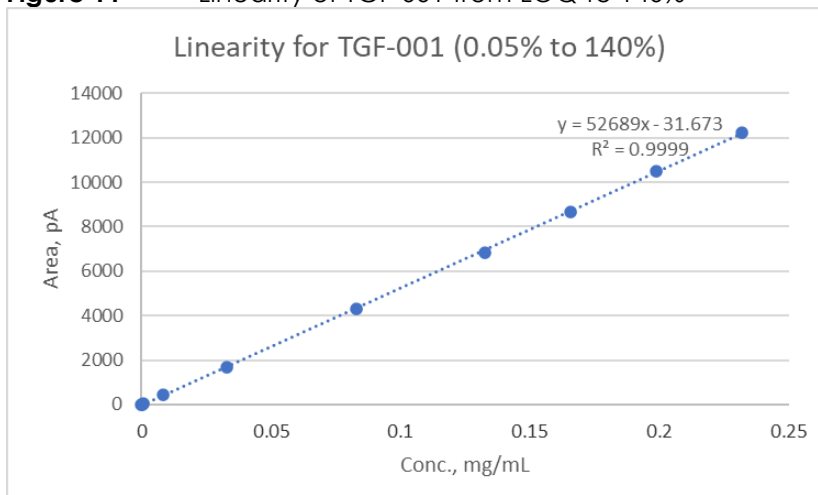
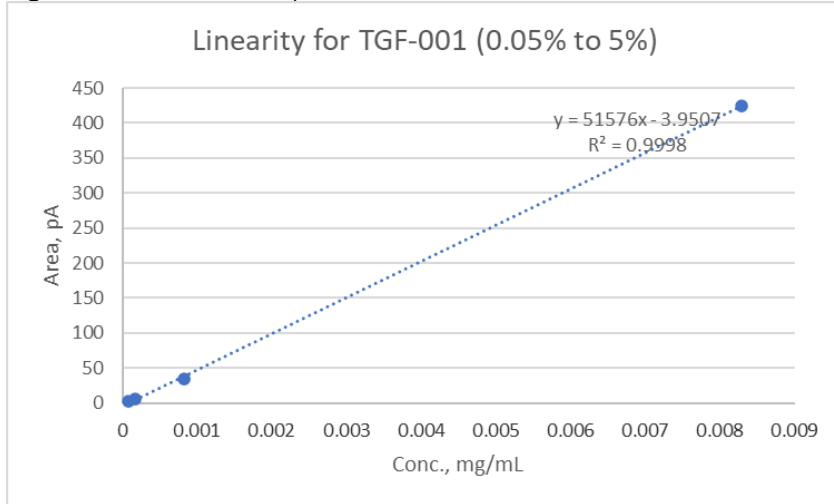
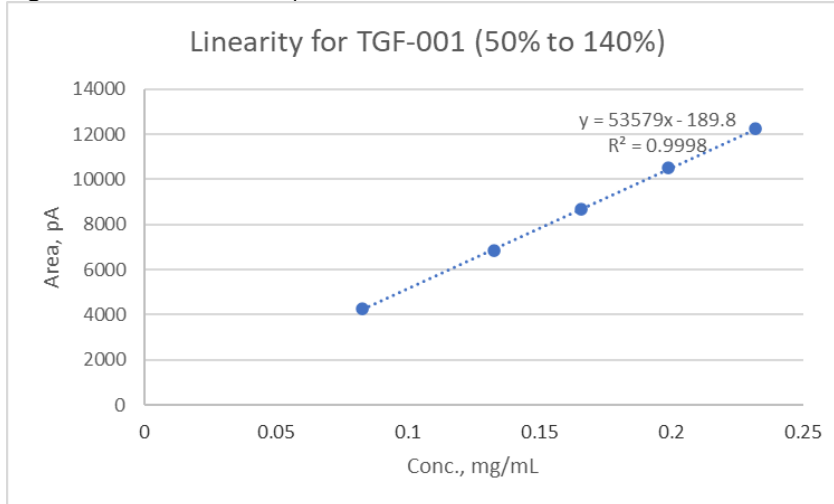


Figure 12 Linearity of TGF-001 from LOQ to 5%**Figure 13** Linearity of TGF-001 from 50% to 140%

3.5. Linearity of DIA and DIN (RRF determination)

Linearity was performed using DIA batch PHTRACKD-394-REF and DIN batch PHTRACKD-389-REF. The correlation coefficient (R^2) was found to be 0.9995 for DIA from LOQ to 0.5% and 0.9993 for DIN from LOQ to 0.5%. All results meet the acceptance criteria.

The Relative Response Factor of DIA relative to TGF-001 was calculated to be '1.16' based on the slope of linearity and could be defined as '**1.00**'.

The Relative Response Factor of DIN relative to TGF-001 was calculated to be '0.91' based on the slope of linearity and could be defined as '**1.00**'.

Table 7 Results for linearity of DIA and DIN

Items	Result	Acceptance criteria
DIA		
R^2	0.9995	≥ 0.995
Slope	45379	report
RRF ⁽¹⁾	1.16	report
DIN		
R^2	0.9993	≥ 0.995
Slope	58071	report
RRF ⁽¹⁾	0.91	report

⁽¹⁾ RRF = slope main compound / slope impurity

Table 8 Concentration and area counts of linearity for DIA and DIN

% of the conc.	Conc. (mg/mL)	Area counts [AU*min]	
		DIA	DIN
0.05%	0.00008	2.3974	4.6605
0.08%	0.00014	4.1521	7.7411
0.1%	0.00017	5.6190	9.3238
0.14%	0.00024	8.5311	13.0105
0.17%	0.00029	10.7812	15.5249
0.3%	0.00051	21.1512	29.4748
0.5%	0.00085	36.6701	49.0316

Figure 14 Linearity of DIA from LOQ to 0.5%


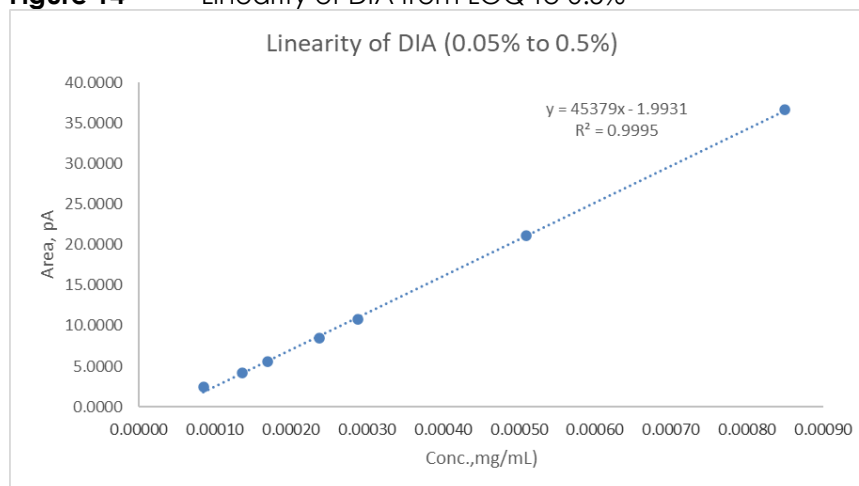


Figure 15 Linearity of DIN from LOQ to 0.5%

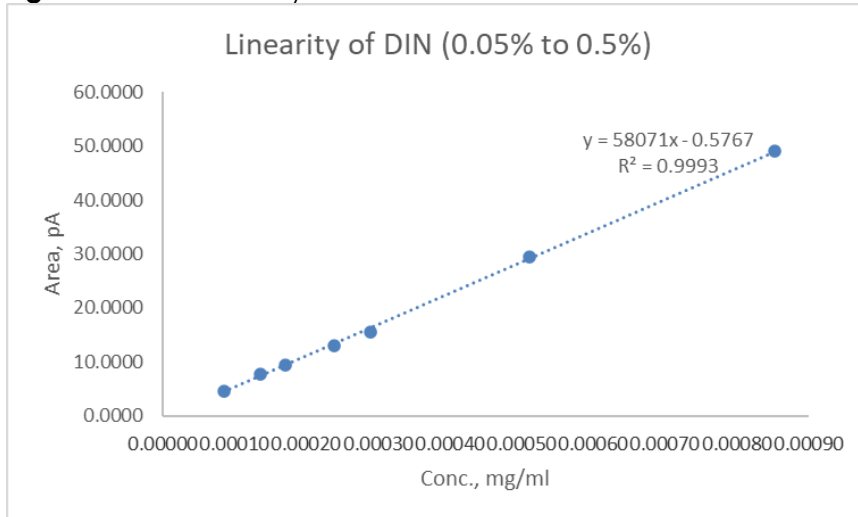

3.6. Accuracy of TGF-001

Accuracy was performed using TGF-001 batch 100144-200502-REF, DIA batch PHTRACKD-394-REF and DIN batch PHTRACKD-389-REF.

The %RSD of TGF-001 peak was determined to be 0.6% at 80% concentration level, 0.2% at 100% concentration level and 1.8% at 120% concentration level. The signal recovery and mean recovery for each concentration level are reported in Table 9, and all results meet the acceptance criteria.

Table 9 Results for accuracy of TGF-001

%RSD for each concentration level	Result	Acceptance criteria
%RSD (80%)	0.6%	≤ 2%
%RSD (100%)	1.7%	≤ 2%
%RSD (120%)	2.0%	≤ 2%
Single recovery	Result	Acceptance criteria
Recovery (80% prep.1)	99%	95% - 105%
Recovery (80% prep.2)	99%	95% - 105%
Recovery (80% prep.3)	98%	95% - 105%
Recovery (100% prep.1)	100%	95% - 105%
Recovery (100% prep.2)	100%	95% - 105%
Recovery (100% prep.3)	97%	95% - 105%
Recovery (120% prep.1)	101%	95% - 105%
Recovery (120% prep.2)	98%	95% - 105%
Recovery (120% prep.3)	98%	95% - 105%
Mean recovery for each concentration level	Result	Acceptance criteria
Recovery (80%)	98%	98% - 102%
Recovery (100%)	99%	98% - 102%
Recovery (120%)	99%	98% - 102%

Table 10 Calculations for accuracy of TGF-001

TGF-001	Weigh in [mg]	Area counts [AU*min]	Response Factor (RF)	Recovery
80% prep.1	11.99	6852.7640	28577.0	99%
80% prep.2	11.96	6821.5891	28518.3	99%
80% prep.3	11.86	6701.9588	28254.5	98%
Average	-	-	28449.9	98%
STDEV	-	-	171.8	-
%RSD	-	-	0.60%	-
100% prep.1	11.99	8660.7480	28893.2	100%
100% prep.2	11.96	8668.1695	28990.5	100%
100% prep.3	11.86	8335.2958	28112.3	97%
Average	-	-	28570.9	99%
STDEV	-	-	481.4	-
%RSD	-	-	1.68%	-
120% prep.1	11.99	10518.3380	29242.0	101%
120% prep.2	11.96	10176.5424	28362.7	98%
120% prep.3	11.86	10029.2460	28187.9	98%
Average	-	-	28597.5	99%
STDEV	-	-	564.9	-
%RSD	-	-	1.98%	-
100% 6 injections (average)	11.99	8663.1630	28901.2944	-

3.7. Repeatability

Repeatability was performed using TGF-001 batch 100144-200502-REF, DIA batch PHTRACKD-394-REF and DIN batch PHTRACKD-389-REF.

The results showed that the %RSD of peak area counts is 0.1% for TGF-001, 0.6% for DIA and 0.9% for DIN. All results meet the acceptance criteria.

Table 11 Results for repeatability for TGF-001, DIA and DIN

100% TGF-001	Area of TGF-001, pA	0.14% DIA in 100% TGF-001	Area of DIA, pA	0.14% DIN in 100% TGF-001	Area of DIN, pA
Inj. 1	8673.119	Inj. 1	13.987	Inj. 1	15.752
Inj. 2	8663.983	Inj. 2	14.085	Inj. 2	15.860
Inj. 3	8665.920	Inj. 3	13.953	Inj. 3	15.506
Inj. 4	8657.123	Inj. 4	14.122	Inj. 4	15.830
Inj. 5	8664.013	Inj. 5	13.972	Inj. 5	15.853
Inj. 6	8654.820	Inj. 6	14.131	Inj. 6	15.802
AVG	8663.1630	AVG	14.042	AVG	15.767
STDV	6.5453	STDV	0.0801	STDV	0.1337
RSD%	0.08%	RSD%	0.57%	RSD%	0.85%
Acceptance criteria	≤ 2%	-	≤ 15%	-	≤ 15%

4. Experimental

4.1. Equipment

- HPLC System: Agilent 1260 Infinity II system
Open Lab CDS-control and integration software or equivalent
- Column: Waters X-Bridge C18, 150*4.6mm, 3.5µm PN: 186003034
- Flow rate: 1.0 mL/min
- Elution: Gradient mode
- Equilibrate time: 10.0 min
- Run time: 40.0 min
- Detection: 278nm
- Injection: 10 µL
- Column temp.: 35°C ± 2°C
- Auto sampler temp.: Room temperature
- Mobile phase:
 - **A:** 0.1% H₃PO₄ in Water
 - **B:** Acetonitrile
- Diluent: 0.1% H₃PO₄ in Water/Acetonitrile (80/20 v/v)
- Needle wash: Water/ Acetonitrile (50:50 v/v)
- Gradient:

Table 12 Gradient Table

Time (min)	% A	% B
0.0	98	2
3.0	98	2
5.0	92	8
25.0	85	15
32.0	15	85
37.0	15	85
37.1	98	2
40.0	98	2

4.2. Equipment and reagents

- Balance: Mettler Toledo XP56
- Acetonitrile: HPLC grade, Merck LiChrosolv, batch K51142730911
- Water: HPLC grade, from Millipore ultra-pure water system
- Phosphoric acid: HPLC grade, Sigmer-Aldrich, batch STBK2887.
- Glassware: 10, 20, 50 mL volumetric flasks, 1L graduated cylinders
- Pipette: 0.2, 1.0 mL, 5.0 mL Pipette

4.3. Solutions

4.3.1. Diluent

0.1% H₃PO₄ in Water/Acetonitrile (80/20 v/v):

Combine 800 mL of water and 200 mL of acetonitrile into a suitable container, and then add 1 mL of Phosphoric acid. Mix well.

4.3.2. Mobile phase

Preparation is described for a volume of 1 liter. Different volumes can be prepared as soon as the solvent ratio is the same.

Mobile phase A (0.1% H₃PO₄ in Water):

In a suitable container, add 1000 mL of water and 1 mL of Phosphoric acid. Mix well.

Mobile phase B (Acetonitrile):

Acetonitrile.

4.3.3. Standard Solutions

➤ Autosampler stability

Accurately weigh approx. 6 mg of TGF-001 reference standard into a 20-mL volumetric flask. Dissolve and dilute to volume with diluent. Mix well.

➤ Specificity/ selectivity and Repeatability

Accurately weigh approx. 6 mg of TGF-001 reference standard into a 20-mL volumetric flask. Dissolve and dilute to volume with diluent. Mix well.

Accurately weigh approx. 6 mg of DIA reference standard into a 20-mL volumetric flask. Dissolve and dilute to volume with diluent. Mix well. Transfer 7 µL of this solution into above 5 mL of TGF-001 solution. Mix well.

Accurately weigh approx. 6 mg of DIN reference standard into a 20-mL volumetric flask. Dissolve and dilute to volume with diluent. Mix well. Transfer 7 µL of this solution into above 5 mL of TGF-001 solution. Mix well.

➤ LOD, LOQ and Linearity

TGF-001 solutions:

Stock Standard Solution (conc.: 0.6 mg/mL, 200%):

Accurately weigh approx. 12 mg of TGF-001 reference standard into a 20-mL volumetric flask. Dissolve and dilute to volume with diluent. Mix well.

140% standard Solution:

Transfer 7 mL stock standard solution into a 10 mL volumetric flask. Fill up to volume with sample diluent and mix well.

120% standard Solution:

Transfer 6 mL stock standard solution into a 10 mL volumetric flask. Fill up to volume with sample diluent and mix well.

100% standard Solution (conc.: 0.3 mg/mL):

Transfer 5 mL stock standard solution into a 10 mL volumetric flask. Fill up to volume with sample diluent and mix well.

80% standard Solution:

Transfer 4 mL stock standard solution into a 10 mL volumetric flask. Fill up to volume with sample diluent and mix well.

50% standard Solution:

Transfer 2.5 mL stock standard solution into a 10 mL volumetric flask. Fill up to volume with sample diluent and mix well.

20% standard Solution:

Transfer 1 mL stock standard solution into a 10 mL volumetric flask. Fill up to volume with sample diluent and mix well.

5% standard Solution:

Transfer 2.5 mL 20% standard solution into a 10 mL volumetric flask. Fill up to volume with sample diluent and mix well.

0.5% standard Solution:

Transfer 1 mL 5% standard solution into a 10 mL volumetric flask. Fill up to volume with sample diluent and mix well.

0.1% standard Solution:

Transfer 1 mL 5% standard solution into a 50 mL volumetric flask. Fill up to volume with sample diluent and mix well.

0.05% standard Solution (LOQ solution):

Transfer 0.5 mL 5% standard solution into a 50 mL volumetric flask. Fill up to volume with sample diluent and mix well.

0.03% standard Solution (LOD solution):

Transfer 0.3 mL 5% standard solution into a 50 mL volumetric flask. Fill up to volume with sample diluent and mix well.

DIA solutions:

Stock Standard Solution (conc.: 0.16 mg/mL, 100%):

Accurately weigh approx. 8 mg of TGF-001 reference standard into a 50-mL volumetric flask. Dissolve and dilute to volume with diluent. Mix well.

0.5% standard Solution:

Transfer 0.5 mL stock standard solution into a 100 mL volumetric flask. Fill up to volume with sample diluent and mix well.

0.3% standard Solution:

Transfer 6 mL 0.5% standard solution into a 10 mL volumetric flask. Fill up to volume with sample diluent and mix well.

0.17% standard Solution:

Transfer 3.4 mL 0.5% standard solution into a 10 mL volumetric flask. Fill up to volume with sample diluent and mix well.

0.14% standard Solution:

Transfer 2.8 mL 0.5% standard solution into a 10 mL volumetric flask. Fill up to volume with sample diluent and mix well.

0.1% standard Solution:

Transfer 2 mL 5% standard solution into a 10 mL volumetric flask. Fill up to volume with sample diluent and mix well.

0.08% standard Solution:

Transfer 1.6 mL 5% standard solution into a 10 mL volumetric flask. Fill up to volume with sample diluent and mix well.

0.05% standard Solution (LOQ solution):

Transfer 0.5 mL 5% standard solution into a 50 mL volumetric flask. Fill up to volume with sample diluent and mix well.

0.03% standard Solution (LOD solution):

Transfer 0.3 mL 5% standard solution into a 50 mL volumetric flask. Fill up to volume with sample diluent and mix well.

DIN solutions:

Same as DIA solutions.

➤ **Accuracy**

Stock Standard Solution (conc.: 0.6 mg/mL, 200%):

Accurately weigh approx. 12 mg of TGF-001 reference standard into a 20-mL volumetric flask. Dissolve and dilute to volume with diluent. Mix well.

120% standard Solution:

Transfer 6 mL stock standard solution into a 10 mL volumetric flask. Fill up to volume with sample diluent and mix well.

100% standard Solution (conc.: 0.3 mg/mL):

Transfer 5 mL stock standard solution into a 10 mL volumetric flask. Fill up to volume with sample diluent and mix well.

80% standard Solution:

Transfer 4 mL stock standard solution into a 10 mL volumetric flask. Fill up to volume with sample diluent and mix well.

Prepare in triplicate.